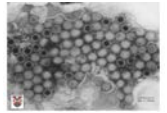
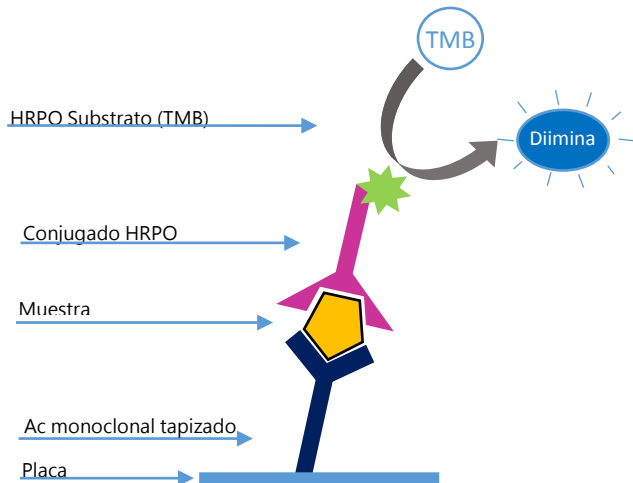
INgezim BTV DAS

R.12.BTV.K2

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
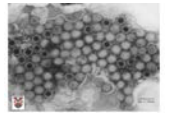
INgezim BTV DAS

R.12.BTV.K2

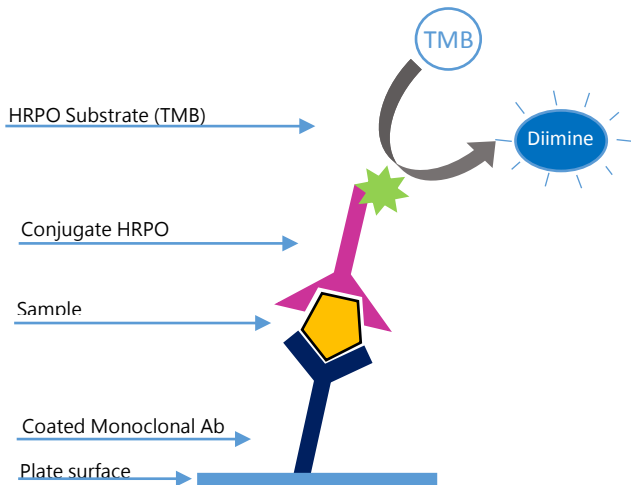INgezim BTV DAS is an enzymatic assay based on a Double Antibody Sandwich ELISA technique, which uses monoclonal antibodies (MAb) specific of Blue Tongue Virus VP7 protein.

TECHNICAL BASIS OF THE KIT

1. Plates are coated with a monoclonal antibody specific of BTV VP7 protein. Samples are added and incubated.
2. If the samples contain BTV, they will bind to the specific monoclonal antibody.
3. After an incubation period, a MAb-HRPO specific to BTV's VP7 is added. It will bind to the protein if there is antigen in the sample.
4. The binding is detected by the development of a colorimetric reaction after the addition of the substrate, specific of HRPO.

APPLICATION

Detection of BTV VP7 antigen in cell cultures and viral load in vaccines.

INTERPRETATION OF THE RESULTS

The amount of VP7 contained in the sample is determined by a Standard Curve obtained for each assay.

VALIDATION

1. Sensitivity of the assay

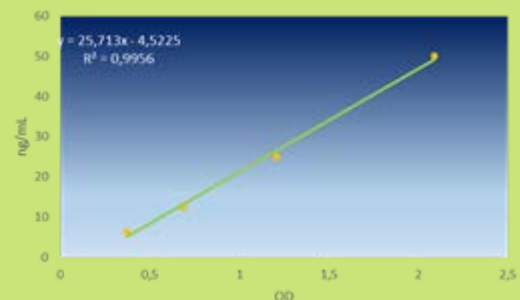
To determine the sensitivity of the assay, different dilutions of a purified VP7 standard were titrated. The results obtained indicated **that the assay is able to detect 3ng of VP7/mL**.

2. Specificity of the assay

To determine the specificity, the assay was tested with African Horse Sickness Virus VP7 protein (virus belonging to the Orbivirus Family). No cross-reaction was observed.

3. Samples quantification

The ELISA DAS can be used to determine precisely the amount of VP7 present in a cell culture sample or in a vaccine, by using a formula which corresponds with a linear regression curve. This curve is obtained with different dilutions of VP7 standard assayed each time.

COMPOSITION OF THE KIT

- Microtitration plate of 96 wells
- VP7 Patron curve
- Vial with Negative Control
- Vial with MAb-HRPO
- Bottle with Washing Solution
- Bottle with diluent
- Bottle with substrate (TMB)
- Bottle with stop solution

PRODUCT MANUFACTURED BY INGENASA

SHELF LIFE: **12 months**
Stored at 2°C-8°C

Ed: 020217